

CARTRIDGE FOR ANALYSIS OF ORGANIC COMPOUNDS IN DRINKING WATER



ENVIRO-CLEAN® METHOD 525.2

ENVIRO-CLEAN® METHOD 525.2

The UCT ENVIRO-CLEAN® Universal cartridge (ECUNI525) provides a new level of performance in solid phase extraction and is specially designed for method 525.2. In addition to its high capture efficiency and broad analyte capture range, use of this fast flow cartridge can significantly increase laboratory sample turnaround time. This brochure contains a new and expanded pesticide list, and updated recovery data including the analyte caffeine.

Product Benefits

- Fast flow rates for rapid analyte capture
- Efficient capture assuring high analyte recovery
- Proprietary bonded C18 with excellent pH stability
- Excellent recovery at all levels of analyte loading
- No lot-to-lot variability
- Excellent reproducibility
- PTFE frits eliminate potential extract contamination
- Cartridge body manufactured from special proprietary polypropylene reducing a potential source of interferences
- Packaged in Mylar to maintain product cleanliness

Product Features

- 83 mL polypropylene cartridge may be used with automated systems
- Can be used with single or multi-station manifolds system
- Each cartridge contains 1500 mg of active sorbent
- Not all C18 bonded silicas can achieve a full range of acceptable recovery using method 525.2. The C18 used in UCT cartridges is specifically designed for 525.2 to enhance the recovery of a wide range of analytes required of the method

Method 525.2 Analytes

The analyte list for this method is comprised of over 120 compounds representative of several classes of pesticides, polynuclear aromatic hydrocarbons, PCBs, phthalates and adipates and other drinking water pollutants. Recovery ranges from 70-130%. Refer to the published method for compound specific MDL's.

The validation data presented herein were determined on independent lots of UCT ENVIRO-CLEAN® Universal Cartridges. MDLs were not determined on all analytes as part of this validation. In addition to the listed method analytes, recovery data for an extended list of analytes is also included.

Table of Compounds Tested Using Method 525.2 and the ENVIRO-CLEAN® Universal Cartridge 525


Analyte	Average 3 Replicates	Std Dev
Acenaphthene	100	0.0
2,4-dinitrotoluene	83	NA
2,6-dinitrotoluene	78	NA
4,4"-DDE	91	4.2
4,4"-DDT	94	3.5
4,4'-DDD	94	4.5
Acenaphthylene	96.0	0.012
Acenaphthene	99.1	0.013
Acetochlor	115	0.01
Alachlor	99	0.007
Aldrin	77	4.4
Ametryn	95	4.6
Anthracene	80	0.0
Atraton	84	17.3
Atrazine	111	0.011
Benzo(a) anthracene	75.4	0.049
Benzo(a)pyrene	105	9.9
Benzo(b)fluoranthene	184.9	0.022
Benzo(k)fluoranthene	95.7	0.029
Benzo[g,h,i]perylene	83.1	0.05
BHC, alpha	108	6.9
BHC, beta	97	3.1
BHC,delta	109	7.9
BHC,gamma	102	11.9
bis- (2-ethylhexyl) adipate	95.1	0.033
bis 2 ethylhexyl phthalate	104	0.029
Bromacil	126	0.012
Butachlor	113	0.005
Butylate	103	4.6
Butylbenzylphthalate	97.1	0.02
Caffeine	90.0	0
Captan	86.9	0.273
Carboxin	103	12.9
Chlordane, alpha	97	4.6
Chlordane, gamma	94	2.5
Chlordane, trans nonachlor	115	11.0
Chlorneb	113	11.0
Chlorobenzilate	118	10.0
Chlorpropham	130	0
Chlorpyrifos (Dursban)	107	5.0
Chlorothalonil	117	12.1
Chrysene	100	0.012
Cyanazine (Bladex)	126	0.008
Cycloate	111	15.0
Dacthal (DCPA) methyl ester	118	13.1
Diazinon	135	0.031
Dibenzo[a,h]anthracene	77.4	0.051
Dichlorvos (DDVP)	127	9.5
Dieldrin	96	6.8
Diethylphthalate	99.1	0.071
Dimethoate	106	0.008
Dimethylphthlate	78.6	0.022
Di-n-butylphthalate	113	0.12

ENVIRO-CLEAN[®] METHOD 525.2

Analyte	Average 3 Replicates	Std Dev
Diphenamid	119	0.008
Disulfoton	92.1	0.01
Disulfoton Sulfone	108	12.5
Endosulfan I	116	11.1
Endosulfan sulfate	114	6.8
Endrin	88	0.0
Endrin Aldehyde	97	3.6
Endrin Ketone	90	3.8
EPTC	102	0.005
Ethion	112	0.005
Ethoprophos	109	5.8
Etridiazole (terrazole)	97	1.2
Fenarimol	70	0.0
Fluoranthene	100	0.018
Fluorene	99.7	0.012
Heptachlor	79	8.2
Heptachlor Epoxide Iso A	116	16.3
Hexachlorobenzene	94	17.4
Hexachlorocyclopentadiene	82	8.4
Hexazinone (Velpar)	105	8.1
Indeno[1,2,3-cd]pyrene	77.4	0.16
Isophorone	91	NA
Lindane	127	4.8
methoxychlor	123	7.6
Methyl Paraxon (Parathion)	115	5.0
Metolachlor	111	0.004
Metribuzin	109	0.005
Mevinphos (phosdrin)	117	12.1
MGK 264	121	5.8
Molinate	114	0.013
Naphthalene	90.3	0.013
Napropamide (Devrinol)	115	2.3
Nonachlor, trans	116	11.1
Norflurazon	133	6.1
PCNB (carbaryl)	91.4	0.021
Pebulate	101	1.7
Pentachlorophenol	80	0.017
Permethrin, cis	124	2.1
Permethrin, trans	123	3.1
Perylene-d12	119	0.0
Phenanthrene	96.9	0.014
Phenanthrene-d10	99	6.6
Prometon	78.6	0.008
Prometryn	110	0.012
Pronamide (propyzamide)	101	1.2
Propachlor	113	15.0
Propazine	105	4.6
Pyrene	94.6	0.022
Simazine	91.4	0.005
Simetryn	93	4.6
Stiufos (tetrachlorvinphos)	126	6.9
Thiobencarb	112	0.008
Tebuthiuron	85	33.5
Terbacil	120	3.5
Terbutryn	103	2.3

Analyte	Average 3 Replicates	Std Dev
Triademefon	98	6.9
Tricyclazole	107	5.0
Trifluralin	82	9.7
Trifluran	83	9.2
Trithion (carbofenothion)	101	0.004
Terbufos	95	7.0
Vernolate	107	1.2
Pyrene	90	7.2
Simazine	94	0.0
Simetryn	93	4.6
Stirofos (tetrachlorvinphos)	126	6.9
Tebuthiuron	85	33.5
Terbacil	120	3.5
Terbutryn	103	2.3
Triademefon	98	6.9
Tricyclazole	107	5.0
Trifluralin	82	9.7
Trifluran	83	9.2
Terbufos	95	7.0
Vernolate	107	1.2

PCB Congeners	Average	Std Dev
2-chlorobiphenyl	93	2.3
2,3-Dichlorobiphenyl	113	15.0
2,4,5-trichlorobiphenyl	97	3.1
2,2,4,4-tetrachlorobiphenyl	98	5.3
2,2,3,4,6-pentachlorobiphenyl	104	2.0
2,2,4,4,5,6-hexachlorobiphenyl	103	3.1
2,2,3,3,4,4,6-heptachlorobiphenyl	85	1.2
Octachlorobiphenyl (BZ#200)	79	1.2

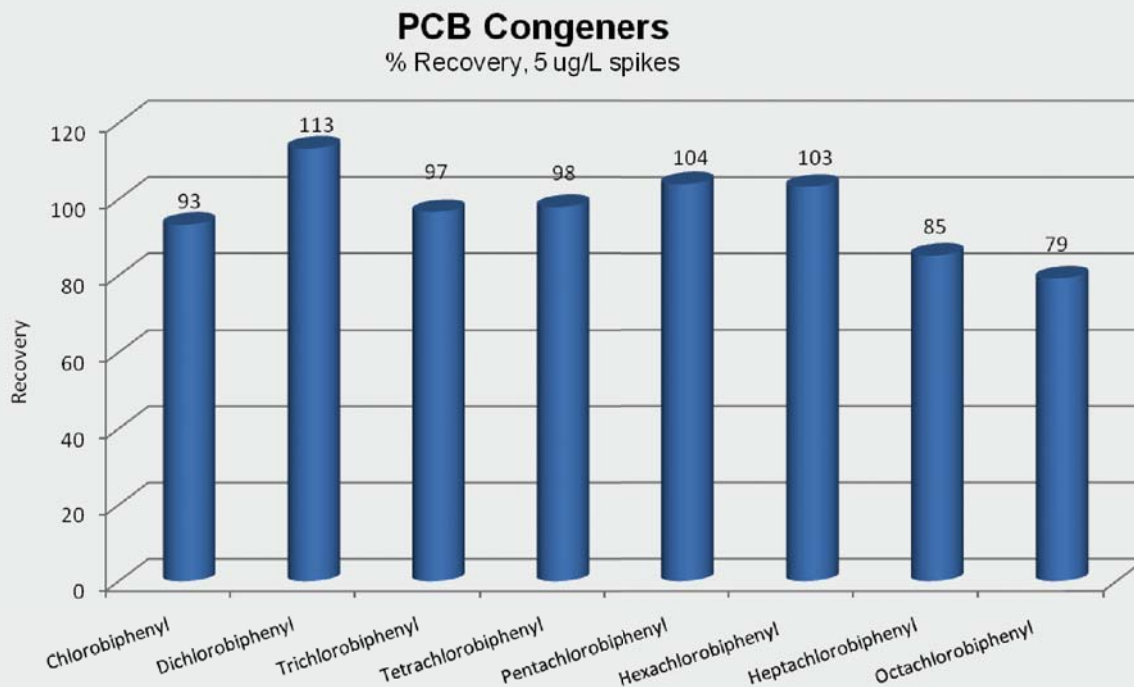

*Data supplied by Rhode Island Department of Health, one replicate. All other data from UCT.

EPA Method 525.2 Revision 2.0

Determination of Organic Compounds in Drinking Water by Liquid-Solid Extraction and Capillary Column Gas Chromatography/Mass Spectrometry

Method Summary

A one liter water sample is adjusted to pH <2 using 6N HCL before passing through an 83 mL ENVIRO-CLEAN® Universal 525 UCT cartridge (ECUNI525). Analytes are eluted from the cartridge with ethyl acetate and methylene chloride. The extract is reduced in volume to 1.0 mL and analyzed by GC/MS.

1. Assemble an all glass filtration apparatus using an 83 mL UCT Universal 525 cartridge.

- a) Use of a vacuum manifold for multiple extractions or automated extractions is acceptable

2. Wash the extraction apparatus and cartridge

- a) Add 10 mL of a 1: 1 mixture of ethyl acetate:methylene chloride (EtAc: MeCl₂) to the reservoir.
- b) Draw a small amount through the cartridge with vacuum
- c) Turn off the vacuum and allow the cartridge to soak for about one minute
- d) Draw the remaining solvent through the cartridge to waste
- e) Allow the cartridge to dry for 3 minutes under full vacuum

3. Condition the cartridge

- a) Add 10 mL of methanol
- b) Draw a small amount through the cartridge
- c) Let soak for about one minute
- d) Draw most of the remaining methanol through the cartridge, leaving 3 to 5 mm of methanol on the surface of the cartridge frit
- e) Immediately add 20 mL of reagent water to the cartridge and draw most of the water through leaving 3 to 5 mm on the top of the cartridge frit

Note: Do not let the cartridge dry out after the addition of water

- f) Add 5 ml of methanol to the water sample and mix well
- g) Add the water sample to the cartridge and under vacuum, filter at a rate of approximately 50 mL per minute
- h) Drain as much water from sample bottle as possible
- i) Dry the cartridge under vacuum for 10 minutes

Note: Exceeding a 10 minute dry time could result in low recoveries. For faster drying, remove the cartridge and tapping the excess moisture from the bottom of the cartridge before continuing vacuum drying.

5. Remove cartridge assembly

- a) Insert a suitable sample tube for eluate collection
- b) Add 10 mL of EtAc to the sample bottle
- c) Rinse the sample bottle thoroughly
- d) Transfer the solvent to the cartridge with a disposable pipette, rinsing sides of filtration reservoir
- e) Draw half of solvent through cartridge then release the vacuum. Allow the remaining solvent to soak the cartridge for about one minute
- f) Draw remainder through under vacuum
- g) Repeat the solvent rinse of the sample bottle and apparatus using 10 mL of 1:1 EtAc:MeCl₂
- h) Using a disposable pipette, rinse down the sides of the cartridge and bottle holder with another 10 mL aliquot of 1: 1 EtAc:MeCl₂
- i) Add the rinse to the cartridge, then draw through

10. Dry the combined eluant

- a) Use granular anhydrous sodium sulfate
- b) Rinse the collection tube and sodium sulfate with two x 3 mL portions of MeCl₂ and place combined solvent in a concentrator tube
- c) Draw through using vacuum
- d) Concentrate the extract to 1 mL under gentle stream of nitrogen (may be warmed gently) being careful not to spatter the contents.

Note: Do not concentrate to <0.5 mL or loss of analytes could occur. Rapid extract concentration could result in loss of low molecular weight analytes

12. Analyze by GC/MS

Revision 2.0, 1995. Method authors: Eichelberger, J. W., Behymer, T. D., Budde, W. L., Munch, J., National Exposure Research Laboratory, Office of Research and Development, U.S. Environmental Protection Agency, Cincinnati, OH 45268

This summary highlights major steps in the 525.2 method. Complete details about the preparation and composition of reagent solutions can be found in method and should be referenced by anyone needing complete details. It is available as a part of Supplement 11 from National Technical Information Service (NTIS), Springfield, VA 22161; publication PB 92 207703. (800) 553-6847 or at www.epa.gov/safewater/methods/methods.html



PRICES AND TERMS

Our prices are subject to change without notice. The price in effect when we receive your order will apply. All prices are in US Dollars and are F.O.B. Lewistown, PA 17044. Terms of payment are net 30 days.

MINIMUM ORDERS

We welcome all orders, therefore, we do not have a minimum order requirement. When ordering, please include your purchase order number, complete "Ship To" and "Bill To" address, catalog number, quantity, and description of product(s). Also include your name and a phone number where you can be reached should we have any questions concerning your order.

SHIPMENTS

Normal processing is within 24 hours after receipt of an order. Unless special shipping requests have been made, our trained staff will send all orders by UPS Ground service. The appropriate shipping charges (freight & insurance costs) will be added to the invoice, unless otherwise instructed by the customer.

SPECIAL PRICING

We offer special pricing for volume purchases and standing orders. These discounts apply to bonded phase extraction column purchases only. Please call a sales representative for more information on special pricing qualifications.

RETURN POLICY

Our Quality Manager will handle all returns. Before returning merchandise, please call to obtain a return authorization number from the quality manager. We will need to know the reason for the return, date of purchase, purchase order number and invoice number in order to issue a return authorization number. Return merchandise must be received before a credit can be issued. Returns will not be accepted after 90 days. A restocking fee of 25% of the price paid, or a minimum of \$25.00 (whichever is greater) will be charged on all returns.

WARRANTY

All products manufactured by UCT are guaranteed against defects in materials and workmanship for a period of 90 days after shipment. UCT will replace any items that prove to be defective during this time period.

The exclusive remedy requires the end user to first advise UCT of the defective product by phone or in writing. Secondly, the defective product must be returned within 30 days after proper approval from our Quality Manager. All returns must indicate the purchase order number, the lot number and the shipping date. UCT's total liability is limited to the replacement cost of UCT products.

This warranty does not apply to damage resulting from misuse.

Placing An Order	Mail	Technical Support
Phone: 215.781.9255 800.541.0559 Fax: 215.785.1226	UCT, LLC 2731 Bartram Rd. Bristol, PA 19007	Phone: 215.781.9255 800.385.3153 Fax: 215.785.1226 Email: info@unitedchem.com Forum: forums.unitedchem.com Web: www.unitedchem.com

