


Barbiturates In Blood, Plasma/Serum, Urine, and Tissue For GC Or GC/MS Confirmations Using: 80 mg Clean Screen® Reduced Solvent Volume Extraction Column

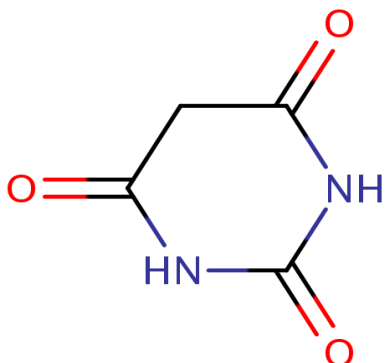

UCT Part Numbers

ZSDAUA08 Clean Screen® RSV - DAU Extraction Column 80 mg 10 mL Without Tips	Or	ZCDAUA08 RSV - DAU w/ Clean Thru® Disposable Tips 80 mg 10 mL
---	----	---

Procedure:

1. Prepare Sample

- To 1 mL of 100 mM phosphate buffer (pH=6.0) add internal standards*. Mix/vortex and add 1 mL of blood, plasma/serum, urine or 1 g (1:4) tissue homogenate.
- Mix/vortex.
- Sample pH should be 5.0 ± 0.5 .
- Adjust pH accordingly with 100 mM monobasic or dibasic sodium phosphate.
- Centrifuge as appropriate.

2. Condition Clean Screen® Extraction Column

- 1 x 0.5 mL CH_3OH .
 - 1 x 0.5 mL D.I. H_2O .
 - 1 x 0.25 mL 100 mM phosphate buffer (pH= 6.0).
- Note:** Aspirate at < 3 inches Hg to prevent sorbent drying.

3. Apply Sample

- Load at 1 to 2 mL/minute.

4. Wash Column

- 1 x 0.5 mL D.I. H_2O ; Aspirate.
- 1 x 0.5 mL 100 mM acetic acid; Aspirate.
- Dry column (5 minutes at > 10 inches Hg).
- 1 x 0.1 mL hexane; Aspirate.

5. Elute Barbiturates

- 1.5 mL hexane/ethyl acetate (50:50).
- Collect at 1-2 mL/minute.

6. Dry Eluate

- Evaporate to dryness at < 40°C.
 - Reconstitute with 100 μL ethyl acetate.
- OPTIONAL DERIVATIZATION
- Add 25 μL of 0.2 M TMPAH****, Reaction occurs in injection port.

7. Quantitate

- Inject 1 to 2 μL onto gas chromatograph.
- For MSD monitor the following ions:



Underivatized				
Compound	Primary Ion**	Secondary	Tertiary	Cerilliant #
Amobarbital	156	141	157	A-020
Butabarbital	156	141	157	B-006
Butalbital	168	167	181	B-024
Hexobarbital*	221	157	236	H-013
Pentobarbital	156	141	197	P-010
Phenobarbital	204	232	117	P-008
Secobarbital	168	167	195	S-002
Thiopental	172	157	173	

Derivatized				
Compound	Primary Ion**	Secondary	Tertiary	Cerilliant #
Butalbital-D5	201	214		B-005
Butalbital	196	195	209	B-024
Amobarbital	169	184	185	A-020
Pentobarbital	169	184	112	P-010
¹³ C ₄ -Secobarbital	200	185		
Secobarbital	196	195	181	S-002
Phenobarbital-D5l	237	151		P-017/P-018
Phenobarbital	232	146	175	P-008

* Suggested internal standard for GC/MS: Hexobarbital or a Deuterated Barbiturate analog.

** Quantitation Ion Target ions in bold

**** Part # STMPAH-0-1, 10, 25, 100

UCT, LLC • 2731 Bartram Road • Bristol, PA 19007 800.385.3153 • 215.781.9255

www.unitedchem.com Email: methods@unitedchem.com

©UCT, LLC 2008 • All rights reserved

