

QuEChERS

Informational Booklet



QuEChERS, the Multiresidue Method of Choice

QuEChERS (pronounced "catchers"), an acronym for Quick, Easy, Cheap, Effective, Rugged and Safe, covers a variety of sample preparation and clean-up techniques for the analysis of multiple pesticide residues in agricultural matrices.

Originally designed for the analysis of fruits and vegetables, **QuEChERS** now includes a wide range of agricultural products. Since its development and publication by scientists at the USDA in 2003, **QuEChERS** has gained significant popularity as the method of choice. It combines several sample preparation steps and extends the range of analytes recovered over older, tedious extraction methods. A driving force in the growth of **QuEChERS** is the emerging need to determine trace amounts of analytes in a high throughput environment.

Matrices include:

- animal products--meat, fish, kidney, chicken, milk, honey
- cereals and grain products
- food products--wines, juices, fruit and vegetables

The expansion of the **Quechers** methodology indicates not only its power for sample extraction and clean-up but also addresses the concern about detecting a vast array of pesticides, herbicides, fungicides, antibiotics, and other compounds throughout the entire food supply.

QuECHERS in its basic form involves three steps:

- 1. liquid micro-extraction
- 2. solid-phase clean-up
- 3. LC/MS/MS or GC/MS analysis

Quechers continues to undergo modifications for improved sample preparation in a broad array of analytes in a vast array of matrices. Due to the large number of Quechers methods now published, Quechers is considered an "approach" rather than a "method." Quechers has now become a generic technique with many modifications, each variation is designed to accomplish one thing—quick sample extraction and clean-up.

Modifications to the original **QuECHERS** method have been introduced to:

- increase sample throughput while reducing costs
- minimize degradation of susceptible compounds (e.g. base and acid labile pesticides)
- expand the range of matrices amenable by this approach





The Three Primary QuEChERS Methods

1) Original QuEChERS Method (by Anastassiades, Lehotay, et al)

- Sodium Chloride is used to reduce polar interferences
- Provides the cleanest extraction because it uses fewer reagents
- Does not use acetic acid which may be problematic in GC/MS analysis
- Uses dispersive clean-up procedures

2) AOAC 2007.01

- Employs 1% acetic acid in acetonitrile and sodium acetate buffer to protect base sensitive analytes from degradation
- A USDA study has demonstrated that this method provides superior recovery for pH sensitive compounds when compared to the other two QuEChERS methods
- The approach uses acetic acid in the extraction step. The acetic acid can overload the PSA sorbent used in the clean-up step making it ineffective and possibly causing GC resolution issues

3) EN 15662

- The European method includes sodium chloride to limit polar interferences and several buffering reagents to preserve base sensitive analytes
- Sodium hydroxide used in the citrus step should be avoided as it can add impurities to the extract as well as damage the sorbent used in the clean-up step

Sample Preparation and Extraction

- Freeze samples to -20°C
- Homogenize with dry ice until a free flowing powder is formed
- The sample is then:
- 1) extracted into solvent
- 2) dispersive or cartridge SPE is used for clean-up

Features and Impact

QuEChERS significantly improves laboratory efficiency and throughput. A batch of 20 extracts can be prepared in less than 60 minutes by a single analyst. This procedure requires only a few milliliters of solvent and is capable of generating recoveries of 90-110% with RSD's < 5% for a wide range of GC and LC amenable compounds.

Extraction and Clean-Up

- · Solvent extraction techniques are designed to remove as much analyte from the base matrix as possible
- Solvent selection is important to minimize co-extracting compounds
- Sample clean-up is necessary to reduce interferences
- Interferences can damage analytical instrumentation and complicate analyte identification and quantification

Extraction Reagents and Their Function

Magnesium sulfate, anhydrous—facilitates solvent partitioning and improves recovery of polar analytes

Acetic acid—used to adjust pH

Acetonitrile—organic solvent providing the best characteristics for extracting the broadest range of pesticides with the least number of co-extractables. Amenable for both LC and GC analysis

Buffers—prevents degradation of pH sensitive analytes by maintaining optimal pH

Sodium Chloride—reduces the amount of polar interferences

Clean-up Reagents and Their Function

Aminopropyl –removes sugars and fatty acids. Serves the same function as PSA, but is less likely to catalyze degradation of base sensitive analytes. Aminopropyl has a lower capacity for clean-up than PSA

ChloroFiltr®— polymeric sorbent for selective removal of chlorophyll from acetonitrile extracts without loss of polar aromatic pesticides

C18—removes long chain fatty compounds, sterols and other non-polar interferences

Graphitized carbon black (GCB)—strong sorbent for removing pigments, polyphenols, and other polar compounds: examples of planar (polar aromatic) pesticides which may be removed: chlorothalonil, coumaphos, hexachlorobenzene, thiabendazole, terbufos, and quintozene

Magnesium sulfate anhydrous—removes water from organic phase

Primary Secondary Amine (PSA)—used in the removal of sugars and fatty acids, organic acids, lipids and some pigments. When used in combination with C18, additional lipids and sterols can be removed

QuEChERS Methods Schematic Flow Chart

Step 1 – Extraction Processes **Buffered QuEChERS Original QuEChERS AOAC QuECHERS Anastassiades and Lehotay 2003 AOAC 2007.01** EN 15662 Add 10 mLs of ACN to 10 g Add 15 mLs of 1% HOAc in ACN to Add 10 mLs of ACN to 10 g homogenized/hydrated sample in 15 ml homogenized/hydrated sample homogenized/hydrated sample in in a 50 mL centrifuge tube a 50 mL centrifuge tube a 50 mL centrifuge tube Add ISTD Add ISTD Add ISTD Shake Shake Shake Add 4 g MgSO₄, 1 g NaCl, 1 g Add 4 g MgSO₄ & 1 g NaCl Shake vigorously for 1 minute Na Citrate 2H, O, 0.5 g Add 6 g MgSO, & 1.5 g NaOAc Na2HCitr-1.5H,O Shake vigorously for 1 minute Centrifuge for 5 minutes Centrifuge at >1500 rcf for 1 minute at 5000 rpm Shake vigorously for 1 minute Centrifuge for 5 minutes at 3000 U/min **Step 2 - Dispersive SPE Clean-Up Processes** Transfer 1 mL aliquot of supernatant Transfer 1 mL aliquot of supernatant to a Transfer 1 mL aliquot of supernatant to a dispersive clean-up tube containdispersive centrifuge tube containing 25 to a micro centrifuge tube containing ing MgSO,, PSA (C18, GCB or Chlomg of PSA and 150 mg MgSO₄, (plus 2.5 150 mg $MgSO_4$ and 50 mg PSA. roFiltr can be added for additional or 7.5 mg of GCB to remove pigments) clean-up) Shake for 1 minute Shake for 30 seconds Shake for 30 seconds Centrifuge for 1 minute at 6000 rpm (5 minutes using GCB) Centrifuge at >1500 rcf for 1 minute Centrifuge for 5 minutes at 3000 U/min. Preserve with toluene for GC/MS or Preserve with 5% formic acid in ACN. Transfer 0.5 mL to vial for GC 6.7mM formic acid for LC/MS/MS or LC analysis Analyze by GC/MS or LC/MS/MS Add TPP surrogate Step 2a - Alternative Cartridge SPE Clean-Up Processes GCB graphitized carbon black Rinse cartridge containing PSA and GCB with 5 mL of acetonitrile MgSO, magnesium sulfate anhydrous **ACN** acetonitrile Transfer an aliquot of the supernatant to the cartridge **HOAC** acetic acid Start collection NaCl sodium chloride

Elute with 6 - 12 mL of 3:1 acetone: toluene

Concentrate for GC/MS or concentrate to dryness and reconstitute in 6.7mM formic acid for LC/MS/MS

Na, Citrate sodium citrate tribasic dihydrate

PSA primary secondary amine

TPP triphenyl phosphate

Na₂HCitr sodium citrate dibasic sesquihydrate

Cartridge or Dispersive SPE (dSPE)

- The original QuEChERS Method uses dSPE clean-up because it's quicker, easier, and less expensive than using traditional SPE cartridges
- With dSPE, the quantity and type of adsorbent can be readily adjusted for differing matrix interferences and various analytes
- · dSPE tubes containing ChloroFiltr® can be used to remove chlorophyll without loss of planar analytes
- PSA and graphitized carbon sorbents are available in 6mL SPE cartridges with PTFE® frits
- · Magnesium sulfate and PSA is available in the Quick QuEChERS cartridge format
- Cartridges provide a better clean-up than dispersive SPE



Polymeric Sorbent

- ChloroFiltr® is a polymeric sorbent available exclusively from UCT. It is designed to replace graphitized carbon black (GCB) for the efficient removal of chlorophyll without loss of planar analytes
- ChloroFiltr® has been tested against hundreds of pesticides and herbicides and has been shown to reduce chlorophyll concentration by greater than 82% without loss of planar analytes.



Comparison of Pesticide Recoveries and RSDs Obtained by dSPE Cleanup of Spinach Sample using ChloroFiltr® and GCB (n=4)				
Pesticide	Chlorof	ChloroFiltr® GCB		GCB
	Recovery%	RSD%	Recovery%	RSD%
Carbendazim	87.1	1.0	71.2	4.0
Thiabendazole	93.2	1.9	55.9	2.6
Pyrimethanil	97.3	1.2	85.0	1.2
Cyprodinil	91.2	0.5	79.3	3.1
Diazinon	104.5	2.3	100.0	0.6
Pyrazophos	92.0	0.9	92.7	1.6
Chlorpyrifos	95.6	2.5	96.3	2.1

The recoveries of carbendazim, thiabendazole, pyrimethanil, and cyprodinil were adversely affected by GCB, especially thiabendazole with a much lower recovery of 55.9% compared to 93.2% obtained by **ChloroFiltr®**. Diazinon, pyrazophos, and chlorpyrifos were less or not affected by GCB due to the non-planar side chains in their structures.



QuEChERS Spinach Extract (acetonitrile) Showing Effectiveness of ChloroFiltr®

Why Use UCT QuEChERS Products?

- Pre-packed products save valuable laboratory time for increased lab throughput
- Best selection of QuEChERS products available
- Cleaner extracts from cleaner products
- Excellent lot to lot reproducibility
- Magnesium sulfate is organic free
- Unique ChloroFiltr® sorbent removes chlorophyll from acetonitrile extracts without loss of planar analytes
- UCT offers sorbents in bulk, dispersive, Quick QuEChERS or traditional cartridge format
- Expert QuEChERS technical support
- · Custom made products are available

Contamination Reduced by UCT Products


A few laboratories assemble their own extraction and clean-up products. QuEChERS sorbents usually become contaminated when exposed to air in the typical laboratory.

A study conducted at a USDA laboratory compared commercially assembled QuEChERS products to those assembled in a USDA lab. Bulk anhydrous magnesium sulfate, PSA, and endcapped C18 sorbents provided by UCT were assembled in a USDA laboratory. These lab preps were compared to UCT assembled products from the same lot of bulk sorbents. The ratio of magnesium sulfate, PSA and C18 was 3:1:1 for this test. Products were evaluated on extracts of milk, honey and soybean and the efficacy of the clean-up was determined by GC/MS analysis. Comparisons of the extracts were made by counting the number of peaks above threshold. Results proved that the UCT assembled products provided superior clean-up compared to the products assembled in the lab. The results were confirmed in three different matrices. The extra peaks observed in the lab prepared products were probably caused by contamination from the lab air. UCT assembled products are prepared under controlled manufacturing conditions so the potential for contamination is eliminated.


These results, along with time and labor savings, prove that QuEChERS products preassembled at UCT are cleaner and more cost effective than products assembled in the lab.

UCT prepared products show a significant reduction in background

Honey Extracted with "In-House" Product

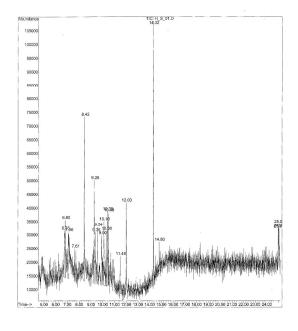

Honey Extract Cleaned with UCT Products

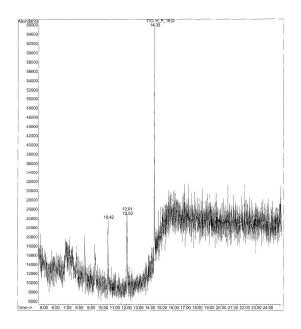

The peaks displayed in the chromatogram on the left show contamination from lab air. The chromatogram on the right shows results from the cleaner UCT prepackaged **Quechers** product.

Studies with soybean and milk products show similar improvement in clean-up when using UCT manufactured vs. laboratory prepared products.

Summary Graph Showing the Total Number of Peaks Seen in GC Chromatograms For Honey, Soybean and Milk



The use of UCT prepared products results in cleaner extracts



I. Recovery Issues

QuEChERS

Tips

- a) Use matrix matched calibration standards for greatest accuracy
- b) Use internal standards
- c) Samples must be at least 80% hydrated for effective extraction
- d) Adding extraction salts directly onto the sample will reduce recovery. Mix sample with solvent first
- e) Buffering is required for base sensitive compounds
- f) Graphitized Carbon Black (GCB) can reduce planar analyte recovery
 - i. Use ChloroFiltr® during extraction to remove chlorophyll
 - ii. Use less GCB
 - iii. Use dual phase (GCB/PSA) cartridge and elute with 3:1 acetone/toluene (product ECPSACB256 is recommended)
- g) For GC analysis, solvent exchanging the final extract into toluene prevents the loss of thermally labile and acetonitrile sensitive pesticides such as chlorothalonil
- h Adding dilute formic acid to the extract after clean-up will prevent degradation of base sensitive compounds while waiting for LC analysis
- i) Do not use **ChloroFiltr®** when extracting mycotoxins. Use endcapped C18 for clean-up.

II. Chromatography Issues

- a) Acetic acid can hinder the clean-up effectiveness of PSA and cause fronting and tailing issues with GC chromatograms. Choose a QuEChERS method that does not use acetic acid
- b) Dispersive SPE may not produce "clean enough" extracts. Use cartridge clean-up to yield a cleaner extract.

Dispersive SPE Clean-Up Guide

- Tube Size Recommendations
 2 mL centrifuge tubes for 1 mL of extract
 15 mL centrifuge tubes for 3+ mL of extract



Matrix	Product Contents	Recommendations Part Number	Product Application & Reference Notes
Pigmented Fruits & Vegetables High pigmentation, some planar analytes	Magnesium sulfate anhydrous Primary Secondary Amine (PSA) Graphitized Carbon Black (GCB) Endcapped C18 (C18)	ECPSAC1856 ECQUEU1115CT ECQUEU32CT ECQUEU515CT ECPSACB256 ECPSACB6 ECPURMPSMC CUMC18CT ECMSC1850CT ECMS12CPSCPSA415CT	F T I W F Z J T U
General Fruits & Vegetables Lightly pigmented	Magnesium sulfate anhydrous Primary Secondary Amine (PSA) Graphitized Carbon Black (GCB) Endcapped C18 (C18)	CUMPS2CT ECMS12CPSA415CT ECMSC1850CT CUMC182CT ECQUEU122CT CUMPSC1875CB2CT CUMPSC18CT ECQUEU1115CT ECQUEU1115CT ECPSACB256	L U T J K, Y K N T
Pigmented Fruits & Vegetables with waxes/lipids	Primary Secondary Amine (PSA) Graphitized Carbon Black (GCB) Endcapped C18 (C18)	ECPSACB6 ECPSAC1856	F F
High Lipid Content (animal products, oils and nuts)	Magnesium sulfate anhydrous Primary Secondary Amine (PSA) Endcapped C18 (C18) Graphitized Carbon Black (GCB)	CUMPS15C18CT CUMPSC1875CB2CT ECMSC1850CT ECPSAC1856 CUMPSC18CT ECQUEU122CT ECPSACB6 EUSILMSSM26 ECMPSC1815CT EEC18156	A, K B, B1, C, C1 F, X X, N, O, P K F G G1 S
Wine and Berries	Magnesium sulfate anhydrous Primary Secondary Amine (PSA) Graphitized Carbon Black (GCB)	ECMPSCB15CT ECPURMPSMC	M M1, Z
Vegetation with Chlorophyll	Magnesium sulfate anhydrous Primary Secondary Amine (PSA) ChloroFiltr® Endcapped C18 (C18)	CUMPSGG2CT CUMPSGGC182CT	R H
Cereal & Grain Products	Magnesium sulfate anhydrous Primary Secondary Amine (PSA) C18 Endcapped (C18)	CUMPS15C18CT EEC18156	D S



UCT QuEChERS Applications Notes

A	Acrylamide by QuEChERS Extraction with LCMSMS Detection	ECMSSC50CT-MP CUMPS15C18CT
В	Flukicides / Anthelmintics by QuEChERS	ECMSSC50CT-MP ECMSC1850CT
B1	Determination of Anthelmintic Drug Residues in Milk Using Ultra High Performance Liquid Chromatography-Tandem Mass Spectrometry	ECMSSC50CT-MP ECMSC1850CT
С	Antibiotics in Beef or Serum by QuEChERS	ECMSC1850CT
C 1	Streamlined Method for the Determination of More Than 100 Veterinary Drugs in Animal Tissue Using Dispersive-SPE Clean-up and LC- MS/MS Detection	ECMSC1850CT
D	Multiresidue Analysis in Cereal Grains Using Modified QuEChERS Method with UPLC-MS/MS and GC-TOFMS	ECMSSC50CT-MP CUMPS15C18CT
E	Trichothecene Type A & B Analysis in Wheat and Corn Using the QuEChERS Approach	ECMSSC50CT-MP CUMPS2CT
F	Modified QuEChERS Procedure for Analysis of Bisphenol A in Canned Food Products	ECQUEU750CT-MP ECPSAC1856 ECPSACB6
G	Determination of Polycyclic Aromatic Hydrocarbons (PAHs) in Seafood Using GC/MS	ECQUUS2-MP ECPAHFR50CT EUSILMSSM26
G1	Extraction of Polycyclic Aromatic Hydrocarbons from Fish Using the QuEChERS Approach	ECMSSC-MP ECPAHFR50CT ECMPSC1815CT
Н	ChloroFiltr®: A Novel Sorbent for Chlorophyll Removal using QuEChERS	ECQUUS2-MP CUMPSGGC182CT
1	Extraction of Pesticides from Tomato Using the QuEChERS Approach This method is applicable to all pigmented fruit and vegetables	ECQUEU32CT ECQUEU750CT-MP ECQUEU515CT
J	Determination of Chlorophenoxyacetic Acid and Other Acidic Herbicides Using a QuEChERS Sample Preparation Approach and LC-MS/MS Analysis	ECQUEU750CT-MP CUMC182CT
К	QuEChERS Sample Preparation For The Analysis Of Pesticide Residues In Olives	ECMSSC50CT-MP ECQUEU122CT CUMPSC1875CB2CT
L	QuEChERS Extraction and Clean-Up of Pesticides from Olive Oil	CUMPS2CT
M	QuEChERS Multiresidue Pesticide Method for the Determination of Multiple Pesticides in Wines	ECQUVIN50CT ECMPSCB15CT

UCT QuEChERS Applications Notes

IV.	VI 1	Determination of Pesticides in Red Wine by QuEChERS Extraction, Quick	ECQUUS2-MP
14		QuEChERS Clean-up, and LC/MS/MS Detection	ECPURMPSMC
N	N	Analysis of 136 pesticides in Avocado using a modified QuEChERS method with LC-MS/MS and GC-MS/MS	ECMSSA50CT-MP CUMPSC18CT
C)	Pesticide Residue Analysis in Whole Milk by QuEChERS and LC-MS/MS	ECMSSA50CT-MP CUMPSC18CT
P		Extraction of Pyrethrin and Pyrethroid Pesticides from Fish Using the QuEChERS Approach	EC4MSSA50CT-MP CUMPSC18CT
C	ם	EURL-FV Multiresidue Method Using QuEChERS by GC-MS/MS & LC-MS/MS for Fruits & Vegetables	ECQUEU750CT-MP ECMPS15CT
R	R	Determination of Pesticide Residues in Marijuana and Tea by QuEChERS and LC/MS/MS	ECQUUS2-MP CUMPSGG2CT
S	8	Analysis of Cyromazine in Poultry Feed Using a QuEChERS Approach	ECMSSA50CT-MP EEC18156
Т	г	QuEChERS Pesticide Analysis for Fresh Produce using GCMSMS	ECMSSC50CTFS-MP ECQUEU1115CT ECMSC1850CT ECMAG00D
L	J	QuEChERS-Based LC/MS/MS Method for Multiresidue Pesticide Analysis in Fruits and Vegetables	EC4MSSA50CT-MP ECMS12CPSA415CT
V	/	QuEChERS Analysis of Miticides and Other Agrochemicals in Honey Bees, Wax or Pollen	ECMSSA50CT-MP CUMPSC18CT ECPSACB256 ECMAG00D
V	N	Multi-residue Pesticide Analysis of Botanical Dietary Supplements using SPE Cleanup and GC-MS/MS	ECMSSC50CT-MP ECPSACB256
×	«	Pesticides in Fatty Matrices	ECPSAC1856 CUMPSC18CT
Y	1	Determination of Carbendazim in Orange Juice Using QuEChERS with LC/MS/MS Detection	ECQUEU750CT-MP CUMPSC18CT
Z	2	Determination of Pesticides in Strawberries by QuEChERS Extraction, Quick QuEChERS Clean-up, and GC/MS Detection	ECQUEU750CT-MP ECPURMPSMC

Products List and Use Description

QuEChERS Multi-Packs

Micro Extraction Products—Reagent Pouches 50 mL centrifuge tubes included (50/pk)

Part Number	Contents
EC4MSSA50CT-MP	4000 mg MgSO ₄
	1000 mg Sodium Acetate
ECMSNA50CT-MP	8000 mg MgSO ₄
	3500 mg Sodium Chloride
EUMIV50CT-MP	6000 mg MgSO ₄
	1500 mg Sodium Chloride
	750 mg Disodium Citrate sesquihydrate
	1500 mg Sodium Citrate tribasic dihydrate
ECMSSA50CT-MP	6000 mg MgSO ₄
	1500 mg Sodium Acetate
ECMSSC50CT-MP	4000 mg MgSO ₄
	1000 mg Sodium Chloride
ECMSSC50CTFS-MP	6000 mg MgSO ₄
	1500 mg Sodium Chloride
ECQUVIN50CT-MP	8000 mg MgSO ₄
	2000 mg Sodium Chloride
ECQUEU750CT-MP	4000 mg MgSO ₄
European QuEChERS	1000 mg Sodium Chloride
Method EN 15662	500 mg Sodium Citrate dibasic sesquihydrate
	1000 mg Sodium Citrate tribasic dihydrate
ECMS4MSC550CT-MP	4000 mg MgSO ₄
	FOO many Conditions Chilesials
	500 mg Sodium Chloride

QuEChERS Multi-Packs

Micro-Extraction Products-Reagent Pouches (without tubes)

Part Number	Contents
ECMSSA-MP	6000 mg MgSO ₄
	1500 mg Sodium Acetate
ECMSSC-MP	4000 mg MgSO ₄
	1000 mg Sodium Chloride
ECQUEU7-MP	4000 mg MgSO ₄
	1000 mg Sodium Chloride
	500 mg Sodium Citrate dibasic sesquihydrate
	1000 mg Sodium Citrate tribasic dihydrate
EUMIV-MP	6000 mg MgSO ₄
	1500 mg Sodium Chloride
	750 mg Disodium Citrate sesquihydrate
	1500 mg Sodium Citrate tribasic dihydrate
ECQUUS2-MP	4000 mg MgSO ₄
	2000 mg Sodium Chloride

Extraction Kits

Part Number	Contents
-------------	----------

ECQUEU215CT	15 mL	6000 mg MgSO ₄
50/pk		1500 mg Sodium Acetate
ECQUEU750CT	50 mL	4000 mg MgSO ₄
50/pk		1000 mg Sodium Chloride
European QuEChERS		500 mg Sodium Citrate dibasic sesquihydrate
Method EN 15662		1000 mg Sodium Citrate tribasic dihydrate
ECMSSC50CT	50 mL	4000 mg MgSO ₄
50/pk		1000 mg Sodium Chloride
ECMSSA50CT	50 mL	6000 mg MgSO ₄
50/pk		1500 mg Sodium Acetate
EUMIV50CT	50 mL	6000 mg MgSO ₄
50/pk		1500 mg Sodium Chloride
		750 mg Disodium Citrate sesquihydrate
		1500 mg Sodium Citrate tribasic dihydrate
ECMS4MSC550CT	50 mL	4000 mg MgSO ₄
50/pk		500 mg Sodium Chloride
Designed for		
Acrylamide Extraction		
ECMS4MSC550CT	50 mL	4000 mg MgSO ₄
		500 mg Sodium Chloride
ECQUEU415CT	15 mL	4000 mg MgSO ₄
		1000 mg Sodium Chloride
		500 mg Sodium Citrate dibasic sesquihydrate
		1000 mg Sodium Citrate tribasic dihydrate



ChloroFiltr® Dispersive Products

Part Number	Size	Contents
CUMPSGG2CT 100/pk A dispersive SPE product for removing polar organic acids, some sugars, lipids an chlorophyll. Designed for 1 n aliquot of supernatant		150 mg MgSO4 50 mg PSA 50 mg ChloroFiltr®
CUMPSGGC182CT 100/pk A dispersive SPE product For removing polar organic acids, some sugars, high lipic chlorophyll. Designed for 1 maliquot of supernatant		150 mg MgSO4 50 mg PSA 50 mg endcapped C18 50 mg ChloroFiltr®
ECMPSGG15CT 50/pk Same as CUMPSGG2CT abo except for larger samples. Designed for 3 mL of supernatant		900 mg MgSO4 300 mg PSA 150 mg ChloroFiltr®
ECMSGG15CT 50/pk Designed for 3 mL of supern	15mL atant	900 mg MgSO4 150 mg ChloroFiltr®

Dispersive Products

Part Number	Size	Contents
ECQUEU12CT	2 mL	150 mg MgSO4
100/pk		25 mg PSA
ECQUEU32CT	2 mL	150 mg MgSO4
100/pk		25 mg PSA
		2.5 mg GCB
ECQUEU42CT 100/pk	2 mL	150 mg MgSO4
		25 mg PSA
		7.5 mg GCB
ECQUEU22CT 100/pk	2 mL	150 mg MgSO4
		25 mg PSA
		25 mg endcapped C18

Dispersive Products

Part Number	Size	Contents
CUMPS2CT	2 mL	150 mg MgSO4
100/pk		50 mg PSA
CUMPSCB2CT	2 mL	150 mg MgSO4
100/pk		50 mg PSA
		50 mg GCB
CUMPSC1875CB2CT	2 mL	150 mg MgSO4
100/pk		50 mg PSA
		7.5 mg GCB
		50 mg endcapped C18
CUMPSC18CT	2 mL	150 mg MgSO4
100/pk		50 mg PSA
		50 mg endcapped C18
CUMPS15C18CT	2 mL	150 mg MgSO4
100/pk		150 mg PSA
		50 mg endcapped C18
ECQUEU122CT	2 mL	150 mg MgSO4
100/pk		50 mg PSA
		50 mg endcapped C18
		50 mg GCB
CUMC182CT	2 mL	150 mg MgSO4
100/pk		50 mg endcapped C18
ECMPS15CT	15 mL	900 mg MgSO4
50/pk		150 mg PSA
ECQUEU315CT	15 mL	900 mg MgSO4
50/pk		150 mg PSA
		150 mg endcapped C18
ECQUEU615CT	15 mL	900 mg MgSO4
50/pk		150 mg PSA
		45 mg GCB
ECQUEU515CT	15 mL	900 mg MgSO4
50/pk		150 mg PSA
		15 mg GCB
ECMPSA50CT	50 mL	1200 mg MgSO4
50/pk	45 1	200 mg PSA
ECMPSCB15CT	15 mL	900 mg MgSO4
50/pk		300mg PSA
F084D00404F0T	45	150 mg GCB
ECMPSC1815CT	15 mL	900 mg MgSO4
50/pk		300mg PSA
		150 mg endcapped C18
ECMS12CPSA415CT	15 mL	1200 mg MgSO4
50/pk		400 mg PSA
CUMPSC1815CT2	15 mL	1200 mg MgSO4
50/pk		400 mg PSA
		400 mg endcapped C18

Dispersive Products

Part Number	Size	Contents
ECQUUS215CT	15 mL	1200 mg MgSO₄
50/pk		400 mg PSA
		400 mg GCB
		400 mg endcapped C18
ECQUEU1115CT	15 mL	1200 mg MgSO ₄
50/pk		400 mg PSA
		400 mg GCB
ECMPSA615CT	15 mL	1800 mg MgSO ₄
50/pk		600 mg PSA
ECMNAX15CT	15 mL	900 mg MgSO ₄
50/pk		150 mg Aminopropyl bonded silica
Florida-Modified QuE	ChERS for	
State Program Fruits a	and Vegetables	
ECMSC1850CT	50 mL	1500 mg MgSO₄
50/pk		500 mg endcapped C18

Cartridge Products
Cartridges are available as an alternative to traditional QuEChERS dSPE clean-up 30/pk

Part Number	Size	Contents		
ECPSACB6	6 mL	200 mg Graphitized Carbon Black GCB (top layer)		
		PTFE frit		
		400 mg PSA (bottom layer)		
ECPSACB256	6 mL	(recommended)		
		250 mg Graphitized Carbon Black GCB (top layer)		
		PTFE frit		
		500 mg PSA (bottom layer)		
ECPSACB506	6 mL	500 mg Graphitized Carbon Black GCB (top layer)		
		PTFE frit		
		500 mg PSA (bottom layer)		
ECMSPSACB6	6 mL	750 mg MgSO ₄ (top layer)		
		500 mg PSA		
		250 mg GCB (bottom layer)		
ECNAXCB506	6 mL	500 mg Graphitized Carbon Black GCB (top layer)		
		PTFE frit		
		500 mg Aminopropyl bonded silica (bottom layer)		
ECPSAC1856	6 mL	500 mg endcapped C18 (top layer)		
		PTFE frit		
		500 mg PSA (Bottom layer)		
ECPURMPSMC	Medium	110 mg MgSO4 (top layer)		
	Cartridge	Teflon Frit		
		180 mg PSA (bottom layer)		
EEC18156	6 mL	500 mg endcapped C18		
EUSILMSSM26	6 mL	200 mg sodium sulfate anhydrous (top layer)		
		1000 mg silica gel (bottom layer)		

QuICK QuEChERS

Simple, Fast, Efficient Cartridges for Clean-Up of QuEChERS Extracts

UCT's QuICK QuECHERS push-thru cartridge eliminates the need for shaking and centrifugation of extracts during clean-up, significantly reducing sample processing time


After QuEChERS sample extraction:

- 1. Draw the extract into a disposable syringe
- 2. Push the extract through the cartridge into a sample vial
- 3. Sample is ready for analysis by LC or GC




The UCT QuICK QuEChERS cartridge ECPURMPSMC contains 110 mg anhydrous MgSO₄ and 180 mg of PSA, providing results comparable to traditional QuEChERS but without the need for centrifugation.


Strawberry Supernatant before Clean-up

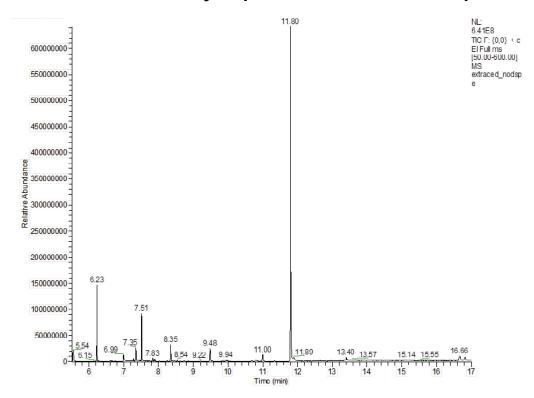

After QuICK Clean-Up

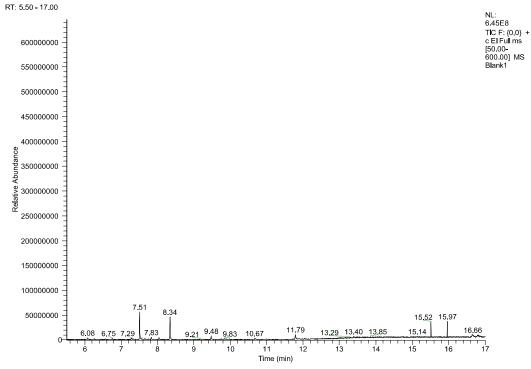

dSPE Recovery Data Comparison vs. QuICK QuEChERS

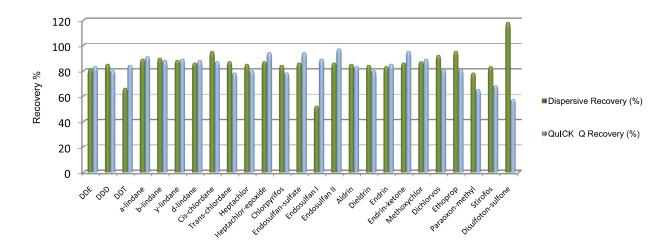



Part Number	Cartridge Size	Contents	
ECPURMPSMC	Medium	110 mg MgSO4 (top layer) PTFE Frit	
		180 mg PSA (bottom layer)	




UCT's new SpinFiltr® is taking the hassle out of dSPE

- Simply discard dSPE chamber containing unwanted matrix and sorbent following centrifugation.
- Recover additional sample volume without worrying about the disruption of centrifugation layers.
- Obtain enhanced extract purification with our built-in 0.2 µm filtration device.

Discover a whole new approach to dSPE!

Product Number	Contents	
ECQUSF14CT	150 mg MgSO₄ 50 mg PSA	
ECQUSF24CT	150 mg MgSO₄ 50 mg PSA 50 mg C18	
ECQUSF34CT	150 mg MgSO₄ 25 mg PSA 2.5 mg GCB	
ECQUSF44CT	150 mg MgSO₄ 50 mg PSA 50 mg C18 7.5 mg GCB	
ECQUSF54CT	150 mg MgSO₄ 50 mg PSA 50 mg C18 50 mg Chlorofiltr®	
ECQUSF64CT	150 mg MgSO ₄ 50 mg C18 50 mg Chlorofiltr®	
ECQUSF74CT	150 mg MgSO₄ 50 mg C18	

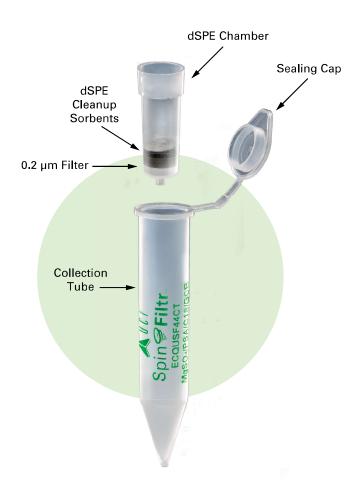

Appendix I

List of possible pesticide analytes that have been shown to yield >90% (or >70 %*) recoveries using the QuEChERS method. GC-amenable pesticides are capitalized; those preferentially analyzed by LC/MS-MS are not capitalized; those that can be analyzed by either technique are underlined**

Pesticide Analytes

acephate*	acetamiprid	Acrinathrin	aldicarb	aldicarb sulfone
aldicarb sulfoxide	Aldrin	azaconazole	azamethiphos	azinphos-methyl
<u>azoxystrobin</u>	Bifenthrin	<u>bitertanol</u>	Bromopropylate	<u>bromuconazole</u>
Bupirimate	<u>buprofezin</u>	butocarboxim	butocarboxim sulfone	butocarboxim sulfoxide
Cadusafos	<u>carbaryl</u>	carbendazim	<u>carbofuran</u>	3-hydroxy-carbofuran
chlorbromuron	(α–, γ–)Chlordane	(α-,β-Chlorfenvinphos	Chlorpropham	Chlorpyrifos
Chlorpyrifos-methyl	Chlorthaldimethyl	Chlorothalonil*	Chlozolinate	clofentezine
Coumaphos	cycloxydim*	(λ -)Cyhalothrin	cymoxanil	Cypermethrin
<u>cyproconazole</u>	<u>cyprodinil</u>	(2,4'-4,4'-)DDE	(2,4'-4,4'-)DDT	Deltamethrin
demeton	demeton-O-sulfoxide	demeton-S-methyl	demeton-S-methyl sulfone	desmedipham
Diazinon	dichlofluanid*	Dichlorobenzophenone	<u>dichlorvos</u>	diclobutrazole
Dicloran	dicrotophos	Dieldrin	<u>Diethofencarb</u>	<u>difenoconazole</u>
Diflufenican	<u>dimethoate</u>	dimethomorph	<u>diniconazole</u>	Diphenyl
Diphenylamine	<u>disulfoton</u>	disulfoton sulfone	diuron	<u>dmsa</u>
dmst	dodemorph	α - Endosulfan	-Endosulfan	Endosulfan sulfate
EPN	<u>epoxiconazole</u>	Esfenvalerate	etaconazole	ethiofencarb sulfone
ethiofencarb sulfoxide	Ethion	ethirimo l	Ethoprophos	<u>etofenprox</u>
Etridiazo l e	Famoxadone	fenamiphos	fenamiphos sulfone	<u>Fenarimol</u>
Fenazaquin	fenbuconazole	fenhexamid*	Fenithrothion	<u>fenoxycarb</u>
Fenpiclonil	Fenpropathrin	Fenpropidine	<u>fenpropimorph</u>	<u>fenpyroximate</u>
enthion	fenthion sulfoxide	Fenvalerate	florasulam*	Flucythrinate I & II
Fludioxonil	flufenacet	Flufenconazole	<u>flusilazole</u>	Flutolanil
Fluvalinate	Fonophos	fosthiazate	Furalaxyl	furathiocarb
<u>furmecyclox</u>	Heptachlor	Heptachlor epoxide	Heptenophos	Hexachlorobenzene
<u>nexaconazole</u>	hexythiazox	imaza l il	imidacloprid	Iprodione
provalicarb	isoprothiolane	isoxathion	kresoxim-methyl	Lindane
linuron	<u>Malathion</u>	<u>malathion oxon</u>	Mecarbam	<u>mephosfolan</u>
Mepronil	Metalaxyl	metconazole	methamidophos*	Methidathion
<u>methiocarb</u>	methiocarb sulfone*	methiocarb sulfoxide	methom yl	methomyl-oxime
metobromuron	metoxuron	Mepanipyrim	Mevinphos	monocrotophos
monolinuron	<u>myclobutanil</u>	nuarimo l	Ofurace	<u>omethoate</u>
<u>oxadixyl</u>	oxamyl	oxamyl-oxime	oxydemeton-methyl	paclobutrazole
Parathion	Parathion-methyl	<u>penconazole</u>	<u>pencycuron</u>	cis- Permethrin
trans-Permethrin	phenmedipham	o-Phenylphenol	<u>Phorate</u>	phorate sulfone
Phosalone	Phosmet	Phosmet-oxon	phosphamidon	Phthalimide
oicoxystrobin	Piperonyl butoxide	<u>pirimicarb</u>	pirimicarb-desmethyl	Pirimiphos-methyl
orochloraz	Procymidone	<u>profenofos</u>	Prometryn	Propargite
Propham	<u>propiconazole</u>	propoxur	Propyzamide	Prothiofos
oymetrozine*	Pyrazophos	pyridaben	pyridaphenthion	<u>pyrifenox</u>
oyrimethani <u>l</u>	Pyriproxyfen	Quinalphos	Quinoxyfen	Quintozene
sethoxydim*	spinosad	<u>spiroxamine</u>	<u>tebuconazole</u>	tebufenozide
<u> Tebufenpyrad</u>	<u>tetraconazole</u>	Tetradifon	Tetrahydrophthalimide	Terbufos
Terbufos sulfone	thiabendazole	thiacloprid	thiamethoxam	thiodicarb
hiofanox	thiofanox sulfone	thiofanox sulfoxide	thiometon	thiometon sulfone
thiometon sulfoxide	thiophanate-methyl	Tolclofos-methyl	tolylfluanid*	<u>triadimefon</u>
<u>triadimenol</u>	Triazophos	trichlorfon	tricyclazole	tridemorph
trifloxystrobin	trifluminazole	Trifluralin	<u>Triphenylphosphate</u>	vamidothion
vamidothion sulfone	vamidothion sulfoxide	Vinclozolin		

^{**}from "Quick, Easy, Cheap, Effective, Rugged and Safe (QuEChERS) Approach for Determining Pesticide Residues", Lehotay, Steven J., U.S. Department of Agriculture, Agricultural Research Service, Eastern Regional Research Center; 600 East Mermaid Lane; Wyndmoor, Pennsylvania 19038; USA

Further Information

Additional information useful to the analyst planning QuEChERS analysis may be found in the following websites:

UCT, LLC

www.unitedchem.com/

A commercial database of application methods and product information related to QuEChERS and other aspects of solid-phase extraction

www.quechers.com

The original website dedicated to the QuEChERS Technique

Nutrient Data Laboratory Website

www.nal.usda.gov/fnic/foodcomp/search/

A nutritional database supported by the USDA Agricultural Research Service

European Websites

http://ec.europa.eu/food/plant/protection/pesticides/index_en.htm

An extensive website maintained by the Health and Consumer Protection Directorate General in Brussels

http://www.crl-pesticides.eu/docs/public/home.asp?LabID=100&Lang=EN

The Community Reference Laboratories and the National Reference Laboratories of the National Food Institute in Denmark













